



The Way Ahead.™

The Challenge of ChIP Sample Preparations

Experiences from Affymetrix and
Recommendations for the Experimental Design

Matthias Prucha, PhD
Senior Applications Specialist and European Tiling
Array Specialist, Affymetrix Europe

1st European ChIP-on-chip Scientific Forum




Outline of this Talk

- **The Challenge**
 - **Technically very different ChIP applications**
 - **General ChIP sample preparation workflow**
- ChIP experimental design
- Affymetrix ChIP protocol

1st European ChIP-on-chip Scientific Forum



Identification of Altered Gene Regulation

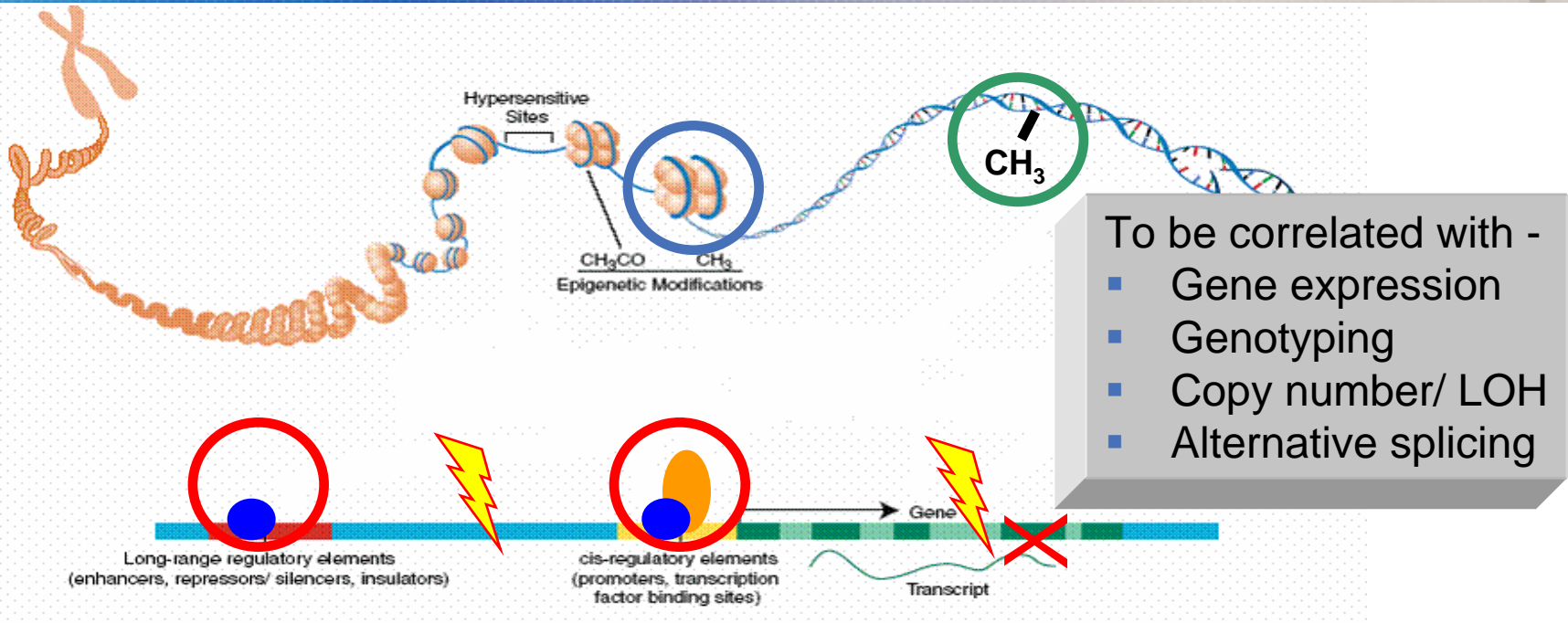

Genome-Wide analysis of transcription regulation using ChIP-on-Chip identifies disease-related modifications at the epigenetic and transcription machinery

- levels:
- **Histone modification**
 - **DNA methylation**
 - **Transcription factor- and other regulatory protein binding activity**
- } Epigenetics

1st European ChIP-on-chip Scientific Forum



(Some) ChIP-chip Publications with Affymetrix Tiling Arrays


- **Cawley S et al., Unbiased Mapping of Transcription Factor Binding Sites along Human Chromosomes 21 and 22 Points to Widespread Regulation of Noncoding RNAs. Cell. 2004 Feb 20;116(4):499-509**
- Bernstein B et al. Genomic Maps and Comparative Analysis of Histone Modification in Human and Mouse. Cell January 28, 2005, Vol 120, 169-181.
- **Carroll J, et al. Chromosome-Wide Mapping of Estrogen Receptor Binding Reveals Long-Range Regulation Requiring the Forkhead Protein FoxA1. Cell, July 12, 2005, Vol. 122, 33-43**
- Schumacher A, et al. Microarray-based DNA methylation profiling: technology and applications. Nucleic Acids research, 2006 Jan., Vol 34 (528-542)
- Bernstein BE, et al. A bivalent chromatin structure marks key developmental genes in embryonic stem cells. Cell. 2006 Apr 21;125(2):315-26
- **Schwartz YB, et al. Genome-wide analysis of Polycomb targets in *Drosophila melanogaster*. Nat Genet. 2006 Jun;38(6):700-5. Epub 2006 May 28. PMID: 16732288**
- Zhang X. et al. Genome-wide High-Resolution Mapping and Functional Analysis of DNA Methylation in *Arabidopsis*. Cell 2006 Sept 22; 126: 1189-1201
- **Carroll J. et al. Genome-wide Analysis of estrogen receptor binding sites. Nat. Genetics Online 2006, Oct 1**
- Hayashi H., et al. High-resolution mapping of DNA methylation in human genome using oligonucleotide tiling array. Hum Genet 2006 Nov, Online

*: In direct relation to Affymetrix ChIP protocol

1st European ChIP-on-chip Scientific Forum



Chromatin Immunoprecipitation Procedure - Optimization Required!

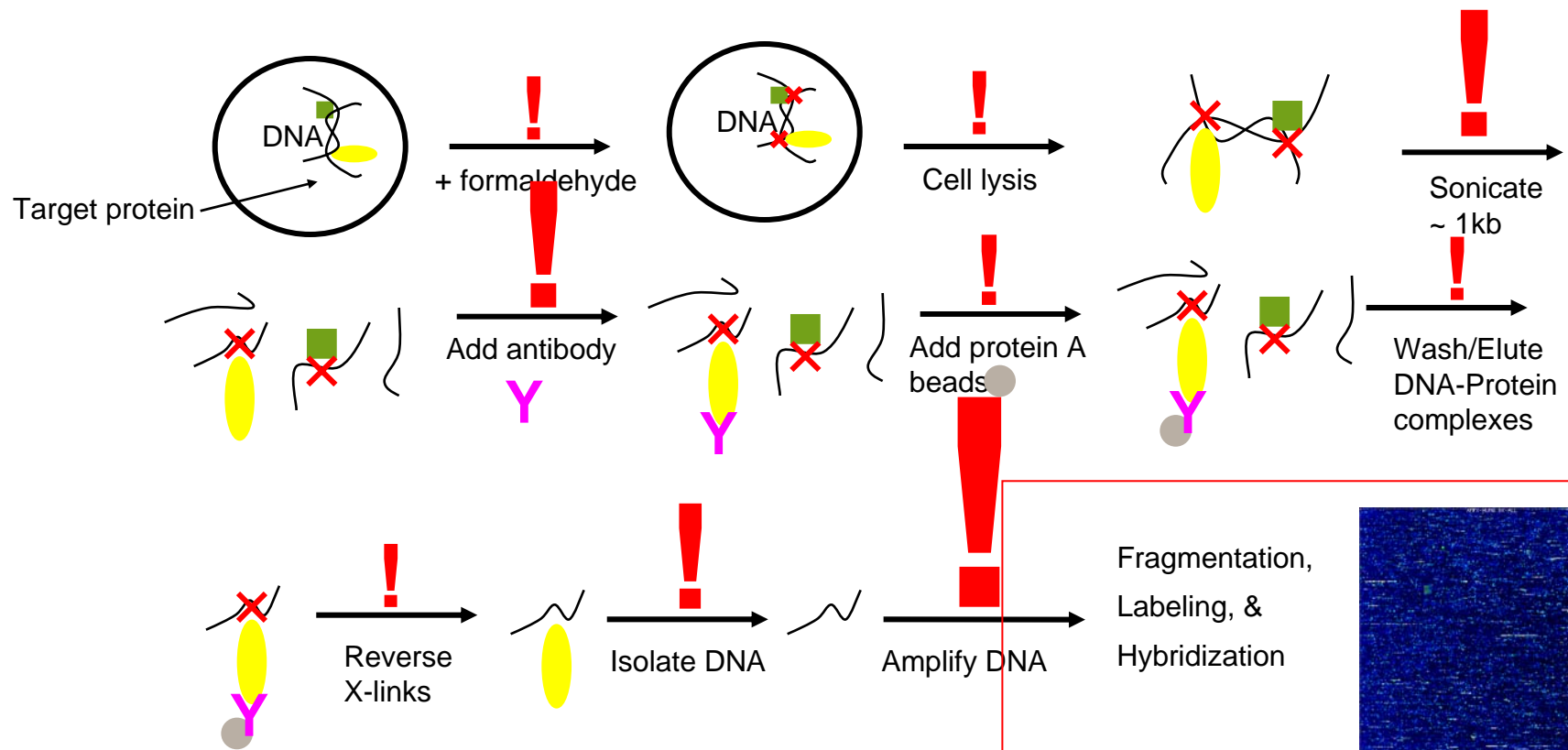

Specific to the Affymetrix technology


1st European ChIP-on-chip Scientific Forum





The Way Ahead.™

The Goal: Hybridization of IP-Enriched DNA

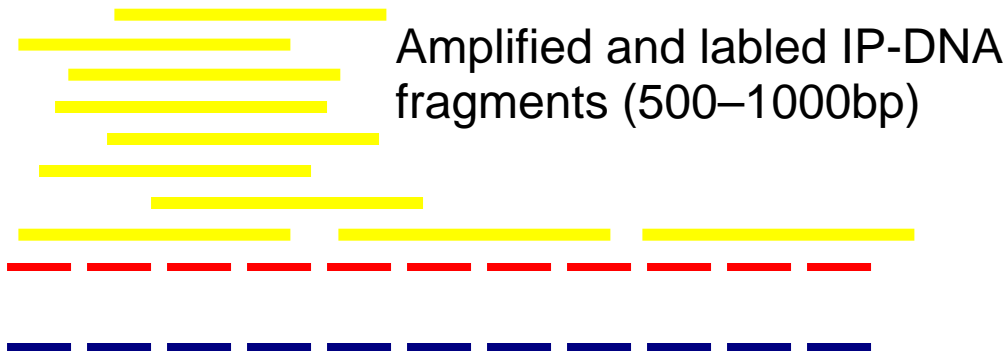


Enriched Sample

PM

MM



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Amplified and labeled non-specific DNA fragments



Control Sample

PM

MM

1st European ChIP-on-chip Scientific Forum



Outline of this Talk

- The Challenge
- **ChIP experimental design**
 - Complexity effects
 - Enrichment analysis
 - Amplification methods
 - Non-specific control
- Affymetrix ChIP protocol

1st European ChIP-on-chip Scientific Forum





The Way Ahead.™

ChIP Challenges are linked!

Complexity

- Genome size
- Frequency of IP sites

Data Analysis

- Normalisation
- Overlap with another specific IP (Poll e.g.)

Antibody

- Affinity
- Specificity
- Stability



Achievable enrichment
(check often, it can get lost)
Enrichment analysis

Hybridization

- Amount per array
- How many cocktails (for array sets)

Amplification

- Which PCR
- Complexity reduction
- Linear amplification (T7)

Non-specific control

- Input
- Mock IP
- Non-specific Antibody

1st European ChIP-on-chip Scientific Forum



Complexity Sources

- **Genome size matters**
 - Human (3200 MB) vs Arabidopsis (157 MB: ~5% of Human)
- **Frequency of binding**
 - Very high
 - DNA-methylation (~19% of entire genome found in Arabidopsis, Zhang et al.)
 - ~Mid to high
 - Histone modifications e.g.
 - ~Low to mid
 - Transcription factors and other regulatory DNA-binding proteins show large differences of binding frequency
- **Antibody-/ purification specificity**
 - Non-specific binding of antibody but also non-specific binding of beads
 - Antibody specificity determines enrichment factor
 - ~95% of the specifically IP'ed DNA is non-specific

1st European ChIP-on-chip Scientific Forum




Complexity Effects

- Amount of labeled DNA for array hybridization
 - Thermodynamics (competition for binding) is correlated with complexity
 - Low complexity (low binding frequency): 0.5 – 10µg per array
 - High complexity: >20µg per array might be useful (see other DNA-applications)
- Selection of amplification method
 - PCR-bias effects (short fragments vs long fragments e.g.) seems to be highly correlated with complexity
 - High complexity: DNA-methylation works in Human and Mouse only with PCR after complexity reduction (Schumacher et al.; GenPathway etc.) or with linear amplification (Hayashi et al.)
 - Low complexity: PCR-methods
- Selection of non-specific IP-control
 - Complexity should be similar between specific IP and unspecific control




Antibody Recommendations

- Only use ChIP-qualified (by Vendor) Ab
- Antibodies vary from batch to batch
 - Use large batches for entire experiment
- Antibody specificity might be unknown
 - Use multiple antibodies for same target
 - Use independent IPs for finding overlaps (bioinformatics approach!)
 - RNA-Polymerase II might be good control for TFs
- Amount of Antibody must be optimized for maximum enrichment (titration)

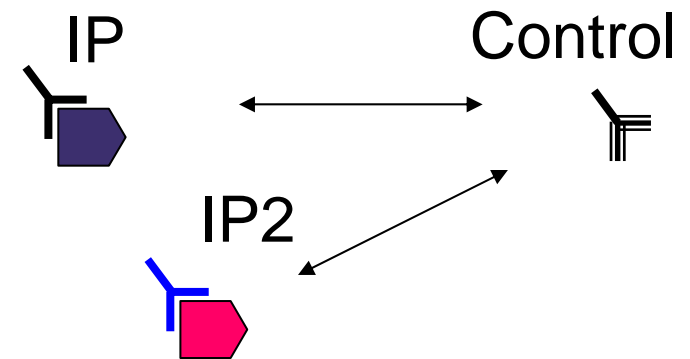

Simplified Workflow – Validation needed

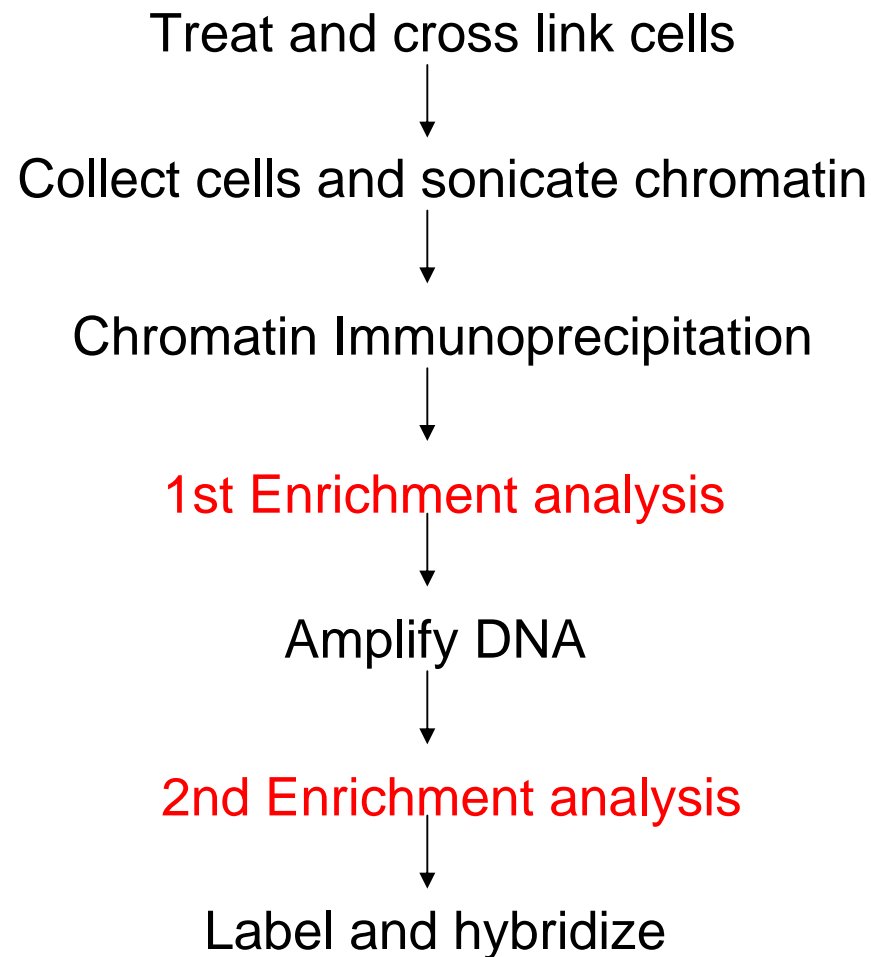


1st European ChIP-on-chip Scientific Forum




Validation – Enrichment Analysis

- Enrichment = sensitivity of ChIP
 - >8-fold is ideal
- PCR of known binding sites
 - Use available hypothesis and literature
 - If unknown binding site... discussed later
- RT PCR versus standard PCR
- Tube to tube variation is guaranteed!
- Need to normalize for differences in DNA concentrations
 - Specific IP vs non-specific control

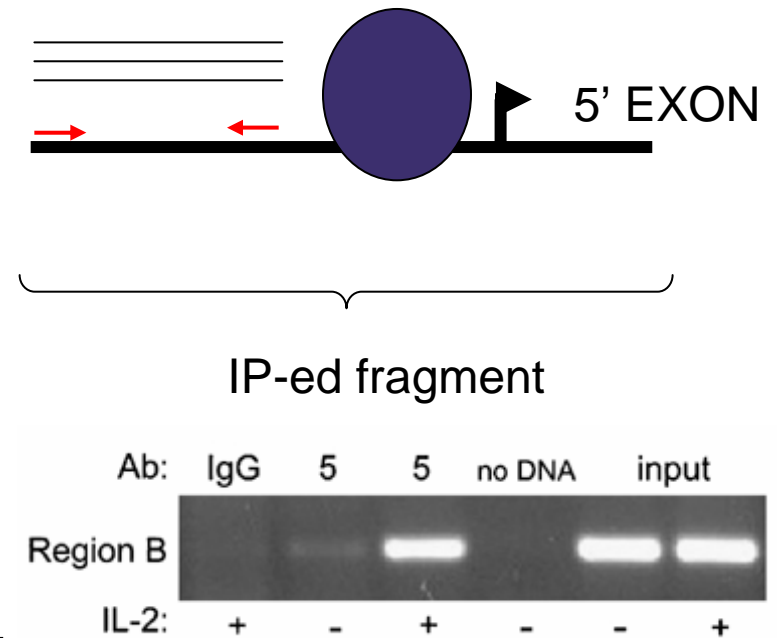

DNA Normalization

- Quantifying DNA is tricky at very low concentrations and large differences between IP- and non-specific control sample
 - Picogreen from Molecular Probes
 - Measures as little 1pg/ul
 - Requires fluorescent plate reader
- Using non-specific region in IP-sample
 - Normalize using primers to a region where no binding occurs
 - Intergenic region, specialized gene (Rhodopsin e.g.)
 - Makes assumption that most ChIP DNA is non-specific

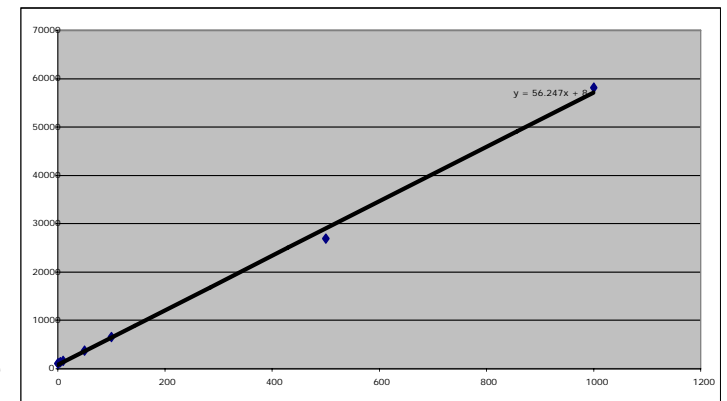

Samples	Fluorescence	pg/ul
ER Chip EtoH	1789	16.66071429
ER Chip E2	1585	13.01785714
Input	62086	1093.392857

1st European ChIP-on-chip Scientific Forum



Unknown Binding Sites

Require a known binding site to monitor through the procedure

- If not:

1. Educated guess. Use expression array data to isolate candidate genes. Make primers to promoter proximal regions
2. Discover binding site: ChIP-cloning approaches
Low success rate, but only need one positive
3. Blind approach: Parallel ChIP of a known positive control (RNA Pol2 etc). If this is successful, then hybridize unknown ChIP sample with one array (one seventh of the genome, chr 21/22, Encodev2)



DNA Amplification – PCR bias?

Many amplification options:



- Random Priming (RP) PCR
 - Ligation Mediated (LM) PCR (used also as complexity red.)
 - Klenow
 - ??? DNA – based assay
- Compatible with Affy ds DNA kit
-
- T7 RP IVT amplification (RNA)
 - Need to be converted to cDNA for hybridization
- Compatible with Affy IVT reagents
- Compatible with Affy ds DNA kit

1st European ChIP-on-chip Scientific Forum



PCR-Saturation affects Enrichment

- Known PCR problems:
 - Bias with fragment length
 - Saturation with too many cycles

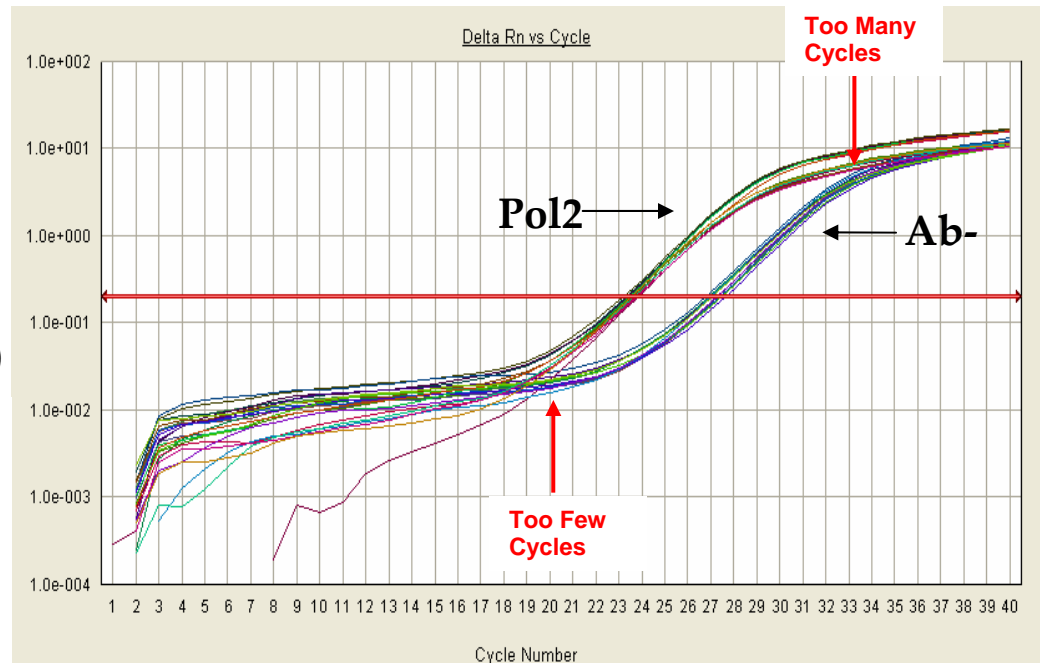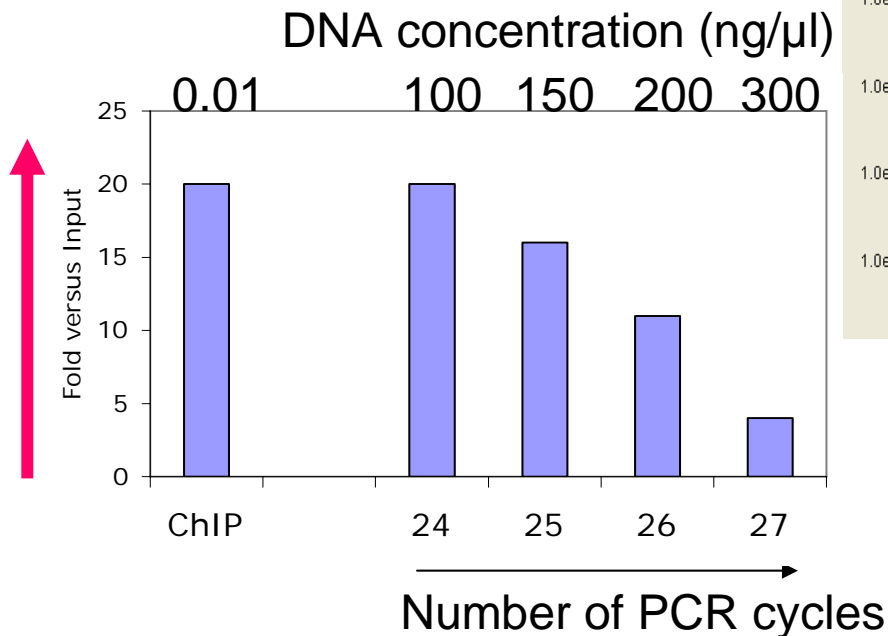

Don't forget to check enrichment after amplification!

1st European ChIP-on-chip Scientific Forum



IP Control Options

- IP experiment (Experimental sample)
 - 10-15ug Ab to IP+ sample (Affymetrix protocol)
- Control experiment
 - Mock IP-DNA (no Ab)
 - Non-specific Ab, e.g. IgG
 - Input-DNA as processed chromatin (X-linked, fragmented)
 - Input-DNA as unprocessed chromatin (genomic DNA)
- Many control options should be tested in parallel.
 - Often Mock IP or using non-specific Ab works best
 - Similar low complexity as specific IP
 - Same concentration range – less error for DNA-normalization
 - Low (or similar) bias during amplification

**Fixed
complexity
(by method)**



**Variable
complexity**



Outline of this Talk

- The Challenge
- ChIP experimental design
- **Affymetrix ChIP protocol**
 - **Sample material, amount of cells**
 - **Cross linking**
 - **DNA fragmentation (shearing)**
 - **Antibody titration, Enrichment analysis**
 - **DNA-Amplification**
 - **Fragmentation and Labeling**
 - **Hybridization; Wash & Stain**

1st European ChIP-on-chip Scientific Forum



Main Message

This protocol is not an
integrated A-Z solution!!!

- The Affymetrix ChIP protocol was developed together with external group at Berkeley University (see Schwartz et al, Mark Biggin – group leader)
- Protocol was validated at Affymetrix by transcription factor binding experiments with Human (SP1, see Cawley et al. 2004) and also Mouse samples
- *Should work for low-mid complexity ChIP applications, higher binding frequencies only in smaller genomes*

1st European ChIP-on-chip Scientific Forum





The Way Ahead.™


Sources of Material

- Suspended cell culture
- Attached cell culture
- Frozen tissues
 - Mouse embryos
 - Drosophila eggs
- Whole blood
- Number of cells
 - Established methods ~ 1×10^8
 - Not practical for many applications
 - Tolerable ~ 1×10^6 cells
 - 1000-10,000 cells using carrier chromatin (O'Neill et al, Nature Genet 2006)
 - Low number of cells leads to loss of enrichment

1st European ChIP-on-chip Scientific Forum



Affy: Cell Preparation / Fixation / Wash

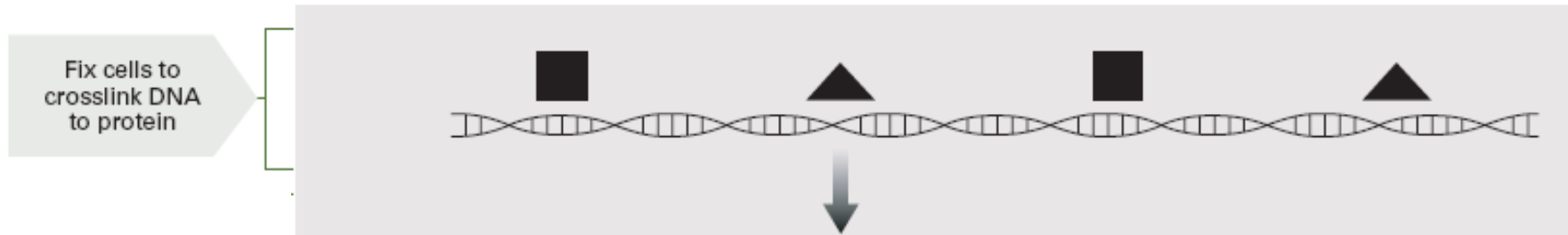


- Cell preparation
 - 5×10^7 /reaction
- Fixation
 - Fix in 1% Formaldehyde 10 min
 - Quench w/ 1/20 vol. of 2.5 M Glycine
- Washing
 - Wash with 1x PBS 2x
 - Wash 3x with Lysis buffer (-80°C Freezing option)

1st European ChIP-on-chip Scientific Forum




DNA Shearing (Fragmentation) Optimization

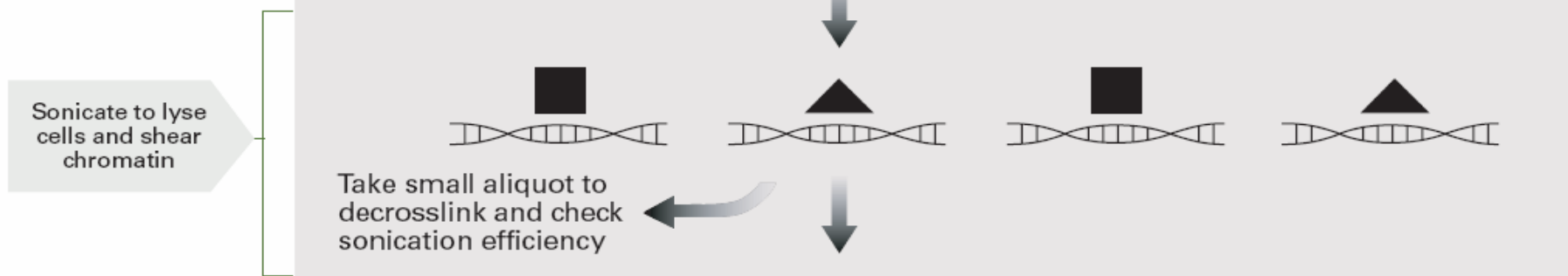

Fragmentation size can also influence PCR (bias effects)

- 1: 2 x 5 sec
- 2: 2 x 10 sec
- 3: 3 x 10 sec
- 4: 3 x 20 sec
- 5: 3 x 30 sec

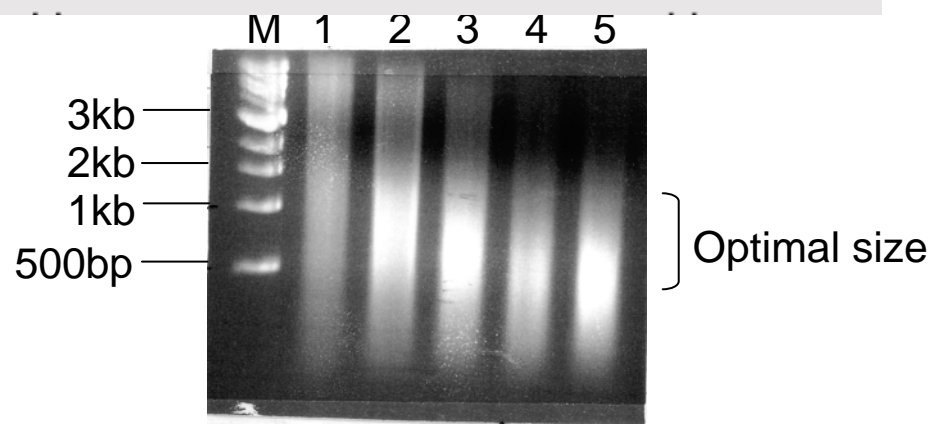


Variables

- Number and duration of pulses
- Sonication needle specs, volume, tube shape and material
- Highly cell type dependent
 - Salt concentrations (Over-fragm. found, lower Ca⁺⁺ and Mg⁺⁺ helped)
 - Additional enzyme (MNase) sometimes needed (Jurkat cells e.g.)



DNA Amplification Step 1 – Linear Amplification

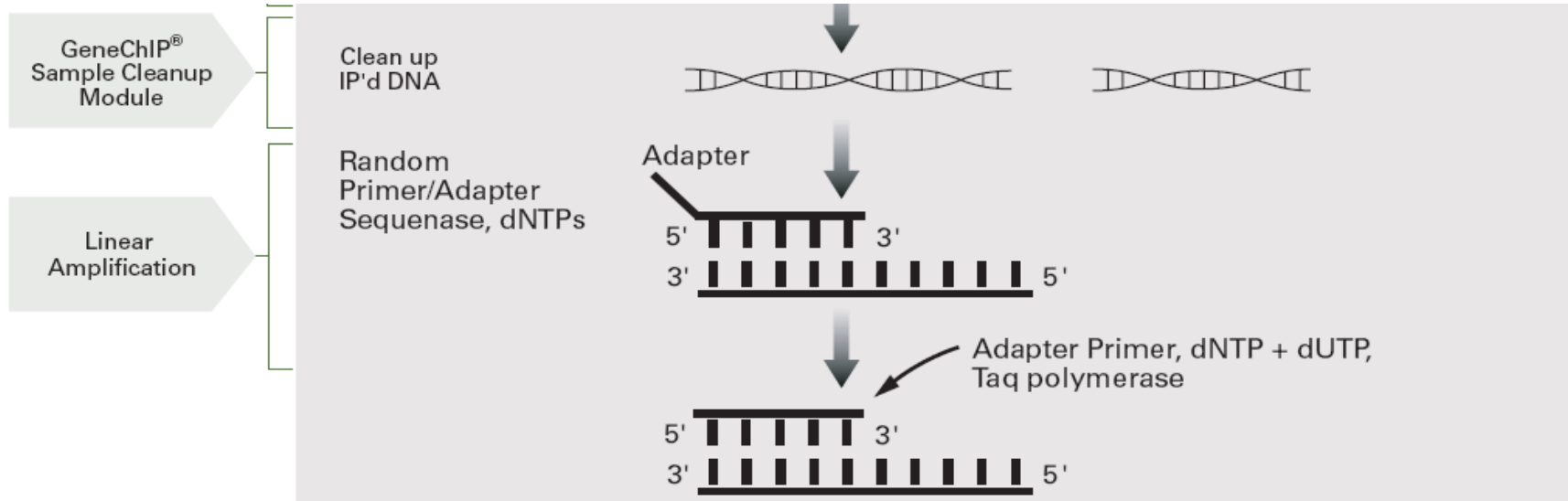

- Tested for low amount of IP-DNA (1-4 ng)
- Substrate inhibition possible
- Temperature gradient set up can be difficult with some Thermocyclers
- Size exclusion columns (Amersham/GEH, two rounds) needed to generate enough yield (remove primers, enrich larger fragments)
 - These columns are storage sensitive


1st European ChIP-on-chip Scientific Forum



Pre-Clearing of Chromatin with Beads

Immunoprecipitate main sample with selected antibody

Couple to Protein-A beads and wash to purify IP'd DNA

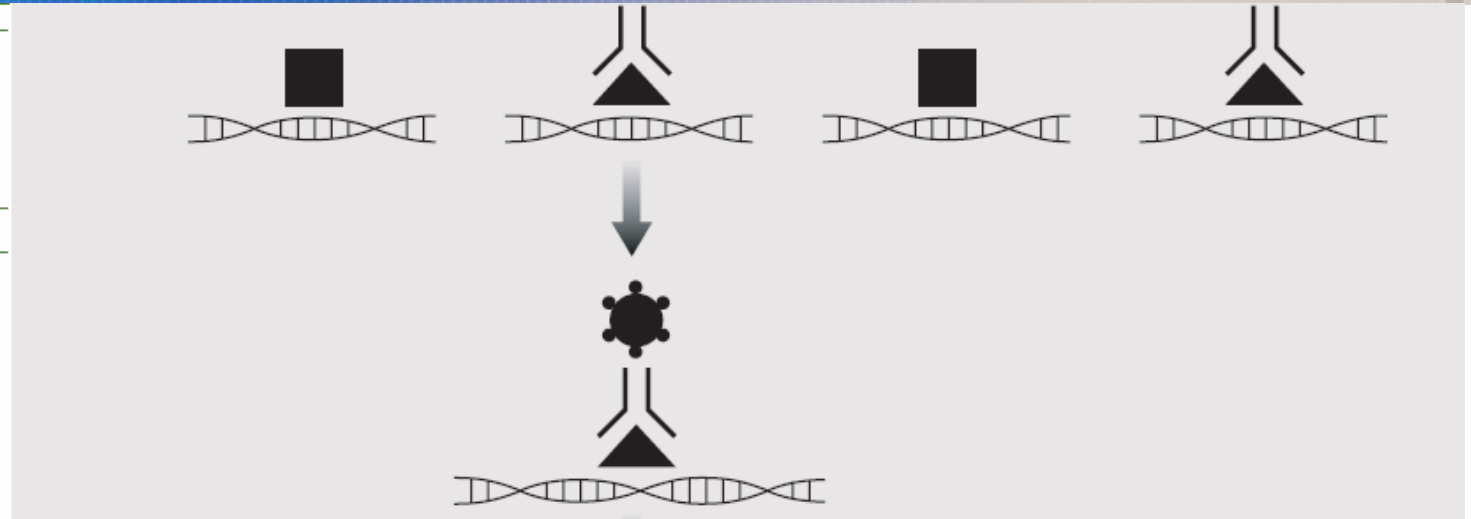


- Part of Affymetrix protocol
- Designed to reduce non-specific interactions of beads with chromatin
 - Some scientists consider it unnecessary
 - Magnetic beads are a good alternative

1st European ChIP-on-chip Scientific Forum




DNA Amplification Step 2: PCR + dUTP Incorporation

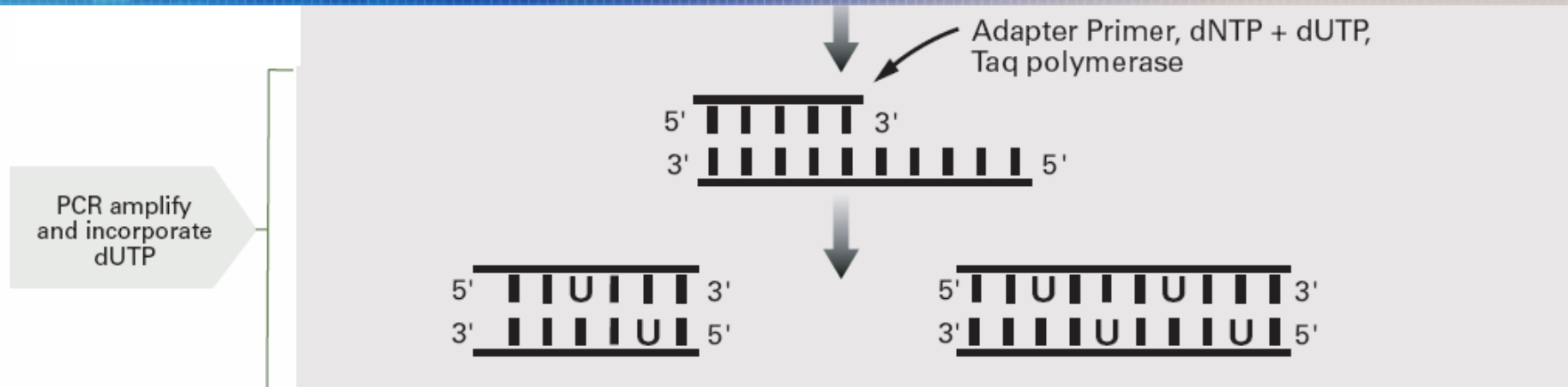

- Problems discussed earlier (complexity-bias; saturation)
- dUTP incorporation prepares for fragmentation reaction
- dUTP can inhibit modern Taq polymerases with side activities (proof reading, often not fully documented)



Fragmentation & Labeling

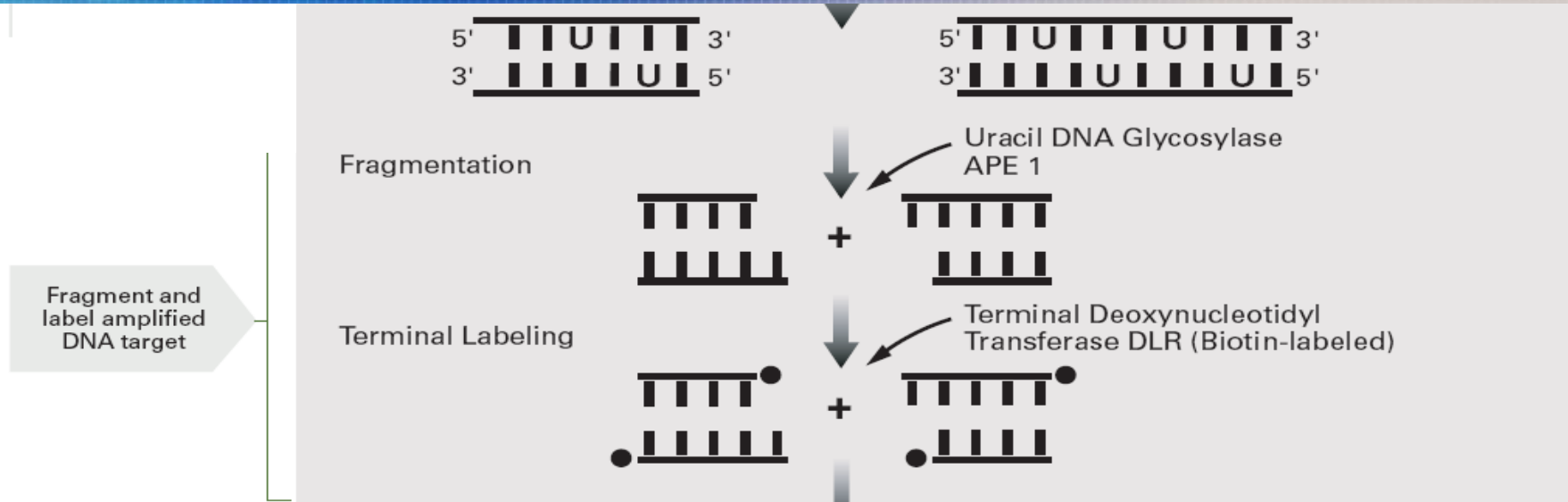


- Fragmentation reaction
 - DsDNA (7-9 ug)
 - Adjust enzymes for DNA amount
 - 37°C 1hour, 93°C 2min
- Labeling reaction
 - Fragmented dsDNA + DLR
 - 37°C 1hour, 70°C 2min

If amplified IP DNA is already available without dUTP used:

- DNA fragmentation & labeling reagents (for 100K SNP-mapping e.g.) can be used as alternative



Hybridization; Wash & Stain

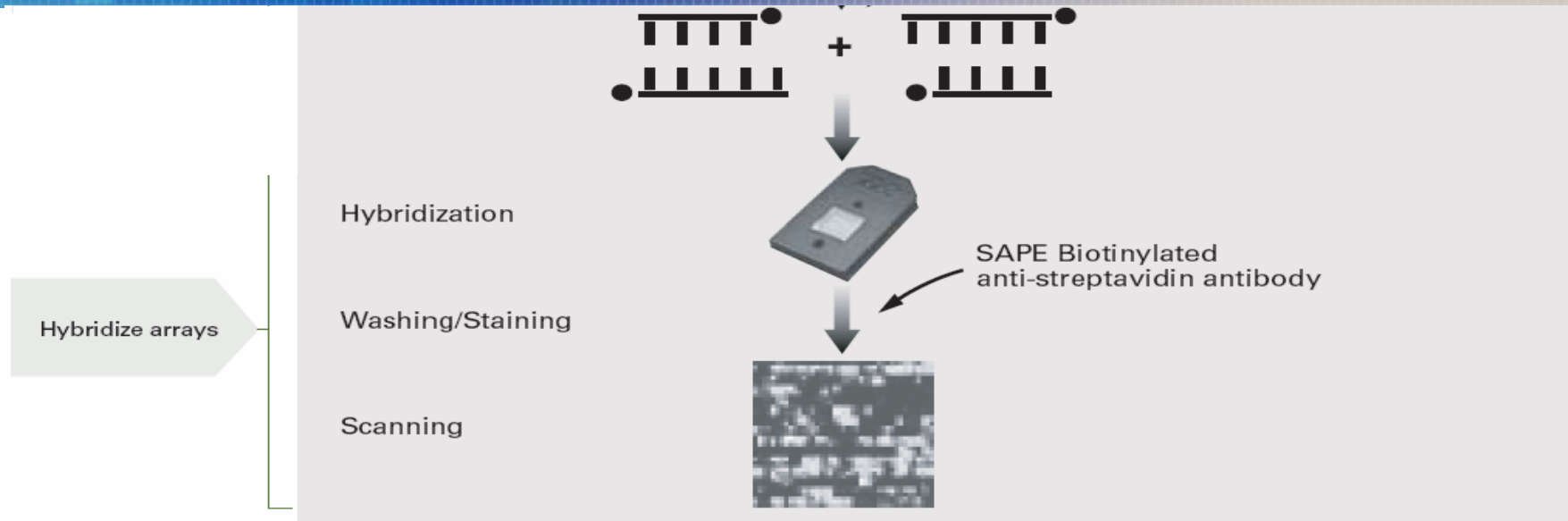

- Amount of fragmented and labeled dsDNA might vary according to complexity and reached enrichment (0.5 – 20 µg)
- Affymetrix Hyb, Wash&Stain kit should be used (w/ additional background blocker)
- 7G GCS-3000 needed ; Specific Fluidics protocols provided

1st European ChIP-on-chip Scientific Forum





The Way Ahead.™

ChIP Assay Optimization (Summary)

- Does intended ChIP application fits to the Affymetrix protocol (complexity...) or do you need to adapt amplification method?
- Sonication conditions of fixed cells
- Antibody titration
 - Antibody affinity and specificity can vary, so the amount of antibody may to be titrated to achieve optimal sample enrichment
- Enrichment analysis needs to be established
 - Repeat after amplification again
- PCR amplification of IP-DNA
 - The optimal number of PCR cycles may require optimization to avoid saturation and ensure that the IP enrichment is maintained
- Adjust amount of amplified dsDNA for fragmentation, labeling and hybridization

1st European ChIP-on-chip Scientific Forum

